

A. Importance Notice

- Please read this manual carefully before use.
- All packaging materials must be disposed of in accordance with local environmental laws & regulations.
- The product shall not be modified or altered in any way. Any modification or alteration will void the warranty.
- Please clean the lid or funnel by following <E. Maintenance> after each use.

B. Unpacking

When unpacking, notice if the following contents are in order. If not, please contact local distributor or email to export@rocker.com.tw for immediate assistance.

Model	PP funnel lid	
Content	PP lid	x1
	Airtight cap	x3
	PTFE Syringe filter	x1
	Instruction manual	x1

C. Installation Example



(a) PP funnel lid




(b) Applied in LF 3



(c) Applied in LF 3a

D. Installation Diagram

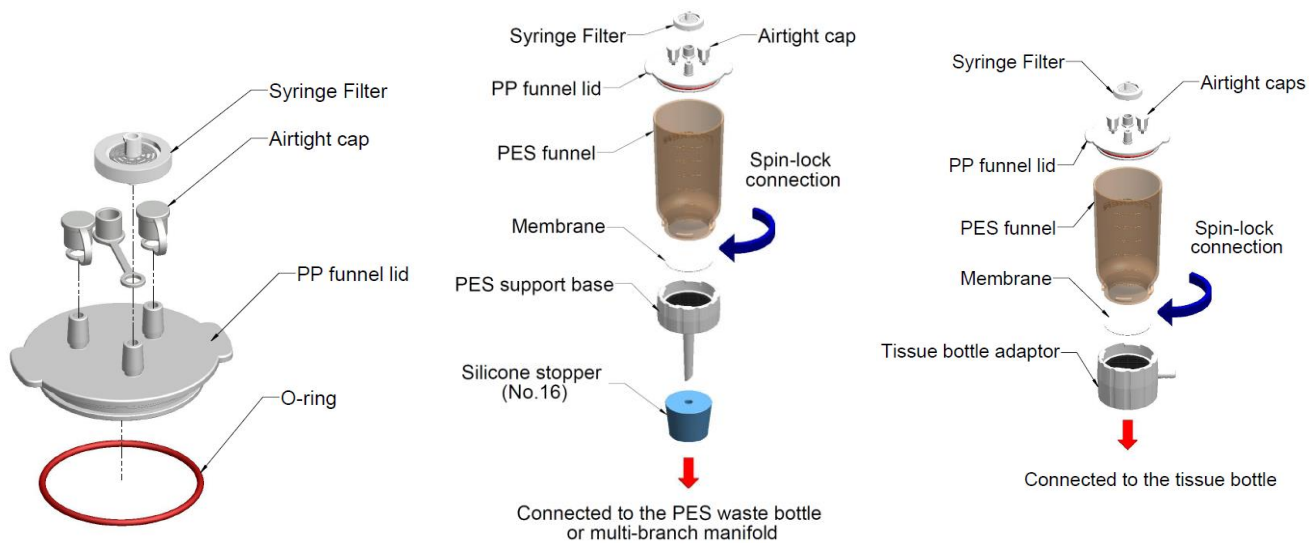

(a) Lid for 300 / 500 ml funnel

(b) LF 3 / 5

(c) LF 3a / 5a

E. Maintenance

1. After each use, clean or sterilize funnel / lid with clean water, 75% alcohol or autoclave.
2. Autoclavable at 121°C for 15 to 20 minutes.

F. Ordering Information

- ◆ 197000-36 PP funnel lid
- ◆ 197000-41 Airtight cap
- ◆ 194225-01 Syringe filter (ø25 mm, 0.2 µm)
- ◆ 197000-44 O-ring for lid kit